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TL-01

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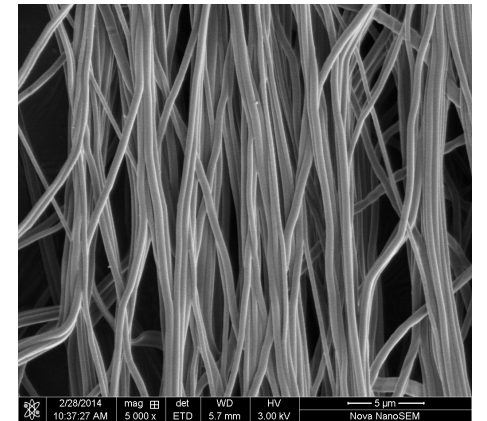
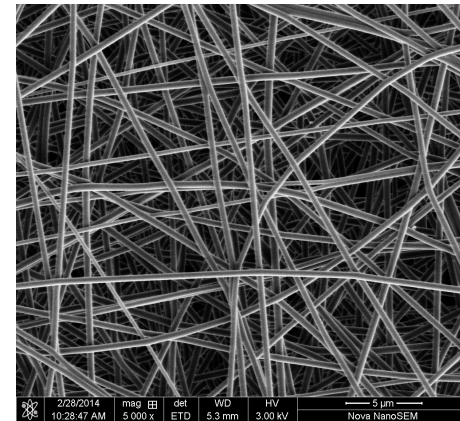
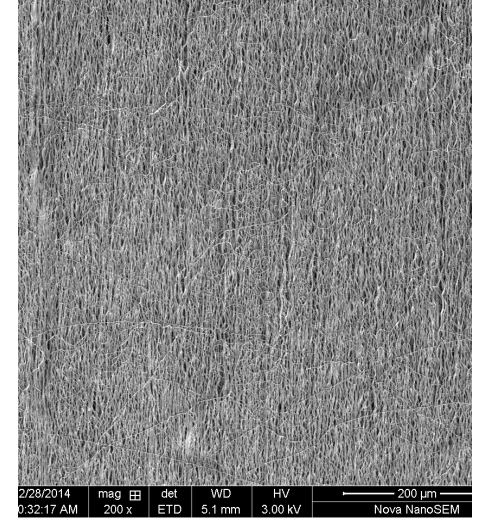
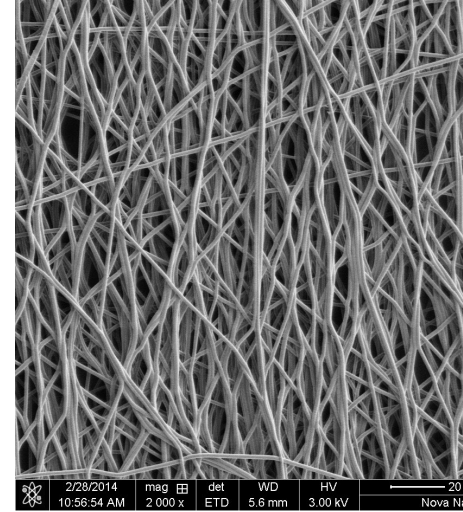
The screenshot displays a control interface with a top status bar showing the date and time '02/23/24 FRI 09:56:52'. Below this, there are fields for 'Now recipe' (0) and 'Now recipe name'. A navigation menu on the left includes 'Constant Mode', 'Variable START', 'Variable STOP', 'View recipe', 'HV', and 'DRUM'. The main control area features 'Homing', 'START', and 'STOP' buttons. A table shows spinning parameters: Spinning time (0 H 0 M 0 S), Rinsing time (0 H 0 M 0 S), and Homing time (0 H 0 M 0 S). On the right, there are 'Heater off' and 'Alarm reset' buttons, and a 'Measured' temperature display showing 0.0. Below this, 'Drum start' and 'Drum stop' buttons are present, along with a table for drum parameters: Drum now speed (0 rpm), Drum setting speed (0 rpm), and ACC/DEC time (0.0 s). Further right, 'Humi' (0.0 %) and 'Temp' (0.0 °C) displays are shown. A bottom navigation bar contains buttons for 'View all parameters', 'Set up', 'Pump', 'Spinning Recipe set', 'Recipe file', 'Information record', and 'System parameters'.

Parameter	Value	Unit
Spinning time	0	H
Spinning time	0	M
Spinning time	0	S
Rinsing time	0	H
Rinsing time	0	M
Rinsing time	0	S
Homing time	0	H
Homing time	0	M
Homing time	0	S
Drum now speed	0	rpm
Drum setting speed	0	rpm
ACC/DEC time	0.0	s

Parameter	Value	Unit
Measured Temp	0.0	°C
Set Temp	0.0	°C
Humi	0.0	%

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